



## Water determination in insulating oil

ESSO UNIVOLT 50 and 56, paraffinic uninhibited transformer oils

HYDRANAL™ Laboratory Report L 462

A coulometric determination was conducted because of the low water content. Since purely alcoholic reagents are not suitable for extracting water with multiple injections, we used reagents that contain solubilizers. While very few oil samples clearly dissolved, they did disperse. Using reagents that contain chloroform and/or xylene requires the use of a coulometry cell with diaphragm.

We found 35 ppm of water with direct injection.

We also followed the indirect method with a KF oven for vials. The analyses showed that 80°C was enough to release the water. We chose to precisely weigh in 4 mL oil sample and we found 26 ppm of water.

In this procedure, the carrier gas is fed directly into the sample. We noted that some 25% of the sample had transferred to the coulometry cell by the end of the determination, because the sample was frothing.

The sample was reduced to around 2 mL so that the carrier gas no longer passed through the sample. This suppressed the frothing. As compensation, we set the minimum determination time to 900 seconds. We found 24 ppm of water.

**Honeywell**

HONEYWELL RESEARCH CHEMICALS PORTFOLIO  
Riedel-de Haën™ Burdick & Jackson™ Fluka™

LABCENTER  
**EXACTA + OPTTECH**



### **Procedure for direct coulometric titration in cell with diaphragm:**

Add approx. 100 mL Hydranal-Coulomat Oil or Hydranal-Coulomat A to the anodic compartment of a coulometric cell with a diaphragm and then add Hydranal-Coulomat CG to the cathodic compartment to the same level (usually 5 mL).

The machine is switched on and it titrates automatically to dryness. When the instrument is showing a low, stable drift, samples can be weighed in by difference using a syringe.

Hydranal-Water Standard 1.0 and Hydranal-Water Standard 0.1 PC are suitable for controlling the coulometric cell.

### **Procedure for indirect coulometric titration with KF oven:**

Place 150 mL Hydranal-Coulomat AG-Oven in the anode compartment of a coulometric cell with diaphragm. Fill the cathode compartment to the same level with Hydranal-Coulomat CG. The cell without diaphragm only requires 150 mL of Hydranal-Coulomat AG-Oven.

The machine is switched on and automatically titrates to dryness. When the drift is low and stable, the carrier gas is connected. When the original stable drift value is approximately reached with the carrier gas, the sample can be weighed in precisely by means of differential weighing and heated to 80°C.

The blank value of the method must be determined and be included in the calculations.

Hydranal-Molecular Sieve 0.3 nm or Hydranal-Humidity Absorber is very well suited as a drying agent for the carrier gas.

Hydranal-Water Standard 1.0 and Hydranal-Water Standard 0.1 PC are suitable for controlling the coulometric cell. Hydranal-Water Standard KF Oven 140–160°C is suitable for controlling the KF oven.



**Europe and International**

**Thomas Wendt**

HYDRANAL Center  
of Excellence

Tel: +49-5137 999-353

Fax: +49-5137 999-698

hydranal@honeywell.com

[34868](#) HYDRANAL-Coulomat Oil

[34807](#) HYDRANAL-Coulomat A

[34739](#) HYDRANAL-Coulomat AG-Oven

[34840](#) HYDRANAL-Coulomat CG

WATER STANDARDS

---

[34828](#) HYDRANAL-Water Standard 1.0

[34426](#) HYDRANAL-CRM Water Standard 1.0

[34446](#) HYDRANAL-Water Standard 0.1 PC

[34693](#) HYDRANAL-Water Standard KF Oven  
140-160°C

AUXILIARIES

---



**Europe and International**

**Agnieszka Kossakowska**

HYDRANAL Technical  
Specialist

Tel: +48 512 355 628

hydranal@honeywell.com

[34241](#) HYDRANAL-Molecular Sieve 0.3 nm

[34788](#) HYDRANAL-Humidity Absorber

HYDRANAL Center of Excellence  
Seelze, June 2003 / October 2017



**USA and Canada**

**Doug Clark**

HYDRANAL Technical Center

Tel: 1-800-Hydranal

(1-800-493-7262)

hydranal@honeywell.com

**For more information, please contact:**

**Exacta+Optech Labcenter Spa**

Via Bosco n.21, 41030 San Prospero (MO)

**Phone:** +39 059 808101

**Email:** info@exactaoptech.com

[www.exactaoptech.com](http://www.exactaoptech.com)

**Honeywell Specialty Chemicals Seelze GmbH**

Wunstorferstrasse 40

30926 Seelze, Germany

Tel.: +49 (0)5137-999-353

Fax: +49 (0)5137-999-698

[hydranal-honeywell.com](http://hydranal-honeywell.com)

All statements and information provided herein are believed to be accurate and reliable, but are presented without guarantee, warranty or responsibility of any kind, express or implied. Statements or suggestions concerning possible use of our products are made without representation or warranty that any such use is free of patent infringement, and are not recommendations to infringe any patent. The user should not assume that all safety measures are indicated herein, or that other measures may not be required. User assumes all liability for use of the information and results obtained. WITHOUT LIMITING THE FOREGOING, HONEYWELL DISCLAIMS THE WARRANTY OF MERCHANTABILITY, FITNESS FOR USE AND NON-INFRINGEMENT. © 2018 Honeywell International Inc.



**RESPONSIBLE CARE**<sup>®</sup>  
OUR COMMITMENT TO SUSTAINABILITY

**Honeywell**

LABCENTER  
**EXACTA + OPTECH**